

This product is **for research use only** (not for diagnostic or therapeutic use)

contact: [support@agrisera.com](mailto:support@agrisera.com)

Agrisera AB | Box 57 | SE-91121 Vännäs | Sweden | +46 (0)935 33 000 | [www.agrisera.com](http://www.agrisera.com)

## Product no **AS15 TMB-HRP-1L**

### AgriseraTMB HRP Substrate (1L)

#### Product information

**Quantity** 1L

**Storage** Stable for 3 years at +2-8° C or at room temperature up to six months, and can also be frozen. Storage at room temperature for 1 month results in ca, 2 % activity loss. Protect the substrate from exposure to sunlight.

#### Additional information

One component ready to use reagent. It contains 2.5 mMol L-1 TMB and Hydrogen Peroxide in a buffer at pH 3.6. It also contains non-toxic stabilizers.

#### How to work with this reagent?

- Pour out amount of reagent which is needed for your assay.
- Do not return any unused substrate back to the original bottle.
- Any trace metals, HRP or other contaminant may be reason for the substrate to turn blue.

#### Application information